

NEW! **accu-jet[®] *pro***

The new BRAND accu-jet[®] *pro* pipette controller builds on the comfort and performance features that made the accu-jet[®] one of the most popular pipette controllers in the market.

- **Even More Comfort.** The accu-jet[®] *pro* has been recontoured for better fit for small and large hands. Weight has been reduced by 10%, and is well balanced for low-fatigue pipetting.
- **Even More Control.** The accu-jet[®] *pro* retains the pressure-sensitive buttons to control pipetting speed found on the previous generation, but adds a motor-speed-limiter control thumbwheel for extra assurance with pipettes as small as 0.1mL.
- **Charge Indicator.** Don't worry about your pipette controller failing on you in the middle of your work. An LED indicates when approximately 2 hours of charge time is left. Instrument can be operated while charging.
- **Advanced Electronics.** The accu-jet[®] *pro* features a long-life, environmentally-friendly Nickel-Metal hydride battery, with a special pulsed-mode charger to eliminate "lazy-battery effect." Instrument supplies approximately 8 hours of continuous pipetting (with a 10mL pipette) without recharging! LED indicates low battery.
- **Ergonomic Design.** The accu-jet[®] *pro* weighs only 6.7oz. (190g) and features a smoothly curved shape to eliminate pressure points for operator comfort.
- **Corrosion Protection.** A built-in check valve protects against liquid penetration. The unique active vapor exhaust system vents externally to protect internal components.



The New Standard in Pipette Controllers.



accu-jet® pro pipette controller

User-accessible battery compartment

Easy-access battery compartment makes changes a snap! Two covers are included - one with ridges to allow inverted storage of the accu-jet® pro on the bench.

Low battery indicator

Red LED lights up when there are two hours of pipetting left in the battery.

Hydrophobic filter

Includes 0.2µm filter for tissue and cell culture applications. Optional 3µm filters are available for pipetting to Class A tolerances.

Silicone Adapter (inside)

The silicone adapter with a built-in check valve provides a firm fit for glass and plastic pipettes from 0.1 to 200mL with a suction tube outer diameter of max. 9.2 mm and reduces risk of over-aspiration. Adapter support, silicone adapter and integral membrane filter are autoclavable at 250°F (121°C).

Mode switch

Select either powered delivery for tissue and cell culture applications, or gravity delivery for analytical applications.

Motor speed control thumbwheel

Adjust maximum pipetting speed for enhanced control with small volume pipettes.

Intuitive push-button control


Press gently for smooth, slow aspiration or dispensing, press more for increased speed. Use in conjunction with the motor-speed control thumbwheel for unsurpassed control for small and large volume pipettes.

Charging socket

The charging socket is conveniently located on the bottom of the grip, keeping the cord out of the way when charging and pipetting simultaneously.

Sculpted grip

Contoured grip eliminates pressure points for comfort in small and large hands.

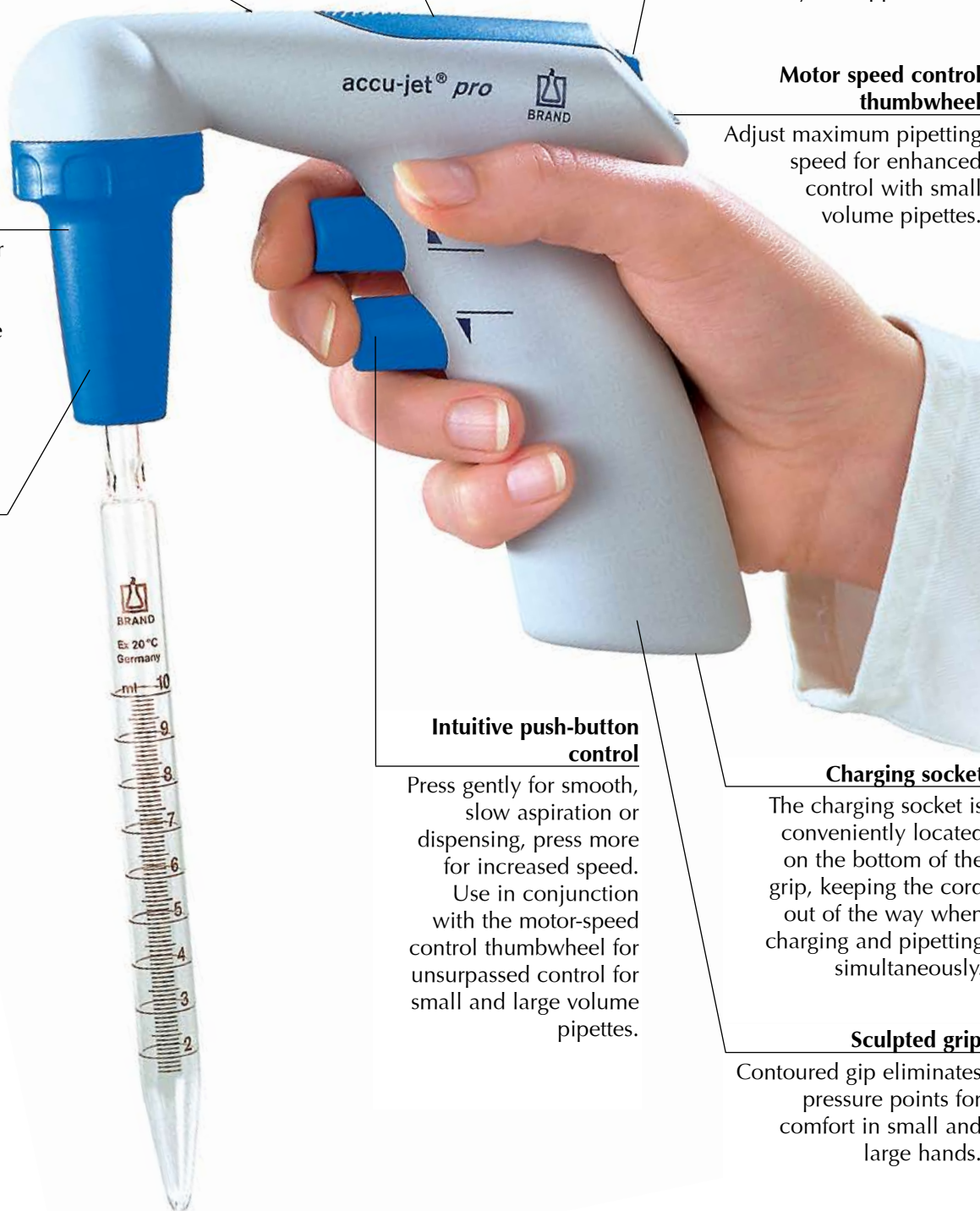

Relaxed Operation

Mission: Control

The accu-jet[®] pro adds a motor speed limiter to the pressure-sensitive buttons to provide accurate, intuitive control for the complete spectrum of pipettes from 0.1mL to 200mL. Two different modes—gravity and powered delivery—greatly expand the range of application.

- **Filling (both models):** Simply adjust the overall pipetting speed using the speed limiter thumbwheel, and use the upper button to bring liquid into the pipette; the powerful internal pump can draw 50mL in less than ten seconds. Aspiration speed is continuously variable by button pressure within the maximum limit set by the thumbwheel. Thumbwheel can be adjusted during use.
- **Gravity-delivery mode:** Free descent of meniscus for volumetric accuracy—ideal for analytical applications.

- **Power-delivery mode:** Continuously adjustable dispensing speed from powered stream to dropwise dispensing including soft blowout—ideal for tissue culture applications.



Recharging

Dual-mode charger provides a quick charging of the environmentally-friendly Ni-MH battery in only four hours. After the full charge is complete, the charger switches automatically to a pulsed mode, to eliminate “lazy battery effect.” The accu-jet[®] pro is fully functional during recharging.



Storage

Ridged battery cover allows inverted storage of the instrument on the benchtop. To save bench space, an included wall support comes with snap tapes for easy cleaning.



NEW! Speed limiter for extra assurance

For extra protection against over-aspiration or excessively powered discharge, set the speed limiter dial at low. For faster or larger volume, set at high. Pipette controller buttons still provide touch-sensitive speed control within the selected speed range.

Technical Data

Items Supplied

The accu-jet® pro pipette controller is supplied with:

- Wall support, gray;
- Nickel-Metal hydride battery;
- Two battery covers; one smooth, one ridged;
- Two spare, sterile 0.2µm membrane filters;
- Charger (120V/60Hz models listed, others available);
- Operating manual;
- One-year warranty.

Operating Limits

Temperature of instrument and reagents must be between 59° and 104°F (15° and 40°C).

Available in colors!

The accu-jet® pro comes in four different color accents for personal preference or color-coding by laboratory or application. Select from dark blue, magenta, green or royal blue.



ORDERING INFORMATION

Description	Cat. No.	2009 List Price
accu-jet® pro Pipette Controller with wall support, spare membrane filters (0.2µm) and charger (120V/60Hz).		
accu-jet® pro Pipette Controller, dark blue	2026330	\$365.00
accu-jet® pro Pipette Controller, magenta	2026331	365.00
accu-jet® pro Pipette Controller, green	2026332	365.00
accu-jet® pro Pipette Controller, royal blue	2026333	365.00
Spare Parts and Accessories		
Wall support, gray	26540	39.00
Silicone adapter with PTFE check valve	26508	23.00
Snap tapes, 12 x 65mm	26542	20.00
Nickel-Metal hydride battery pack	26630	23.00
Charger, 120V/60Hz	26603	81.00
Membrane filter, 0.2µm, blister pack, sterile*, pack of 10	2026539	91.00
Membrane filter, 0.2µm, non-sterile, bag of 10	26535	75.00
Membrane filter for Class A pipetting, 3µm, blister pack, non-sterile, each	26052	11.00
Membrane filter for Class A pipetting, 3µm, non-sterile, pack of 10	26056	61.00

*Sterilized with ethylene oxide (EO), with a guaranteed SAL of 10⁻⁶.